


ezImageX3 Setup Instructions

Whether you are using 1 or 2 cameras it is necessary to complete all the setup steps below to ensure that the program will function as intended. This also ensures that your user settings can be saved.

1. After installing ezImageX3, navigate to the **View** Tab and click **Dual** to enter dual camera mode.




2. With Camera 1 selected, navigate to the **Capture** Tab and click **Edit** to access the edit menu.
3. Identify your camera by device name within the first drop-down list and select that camera.

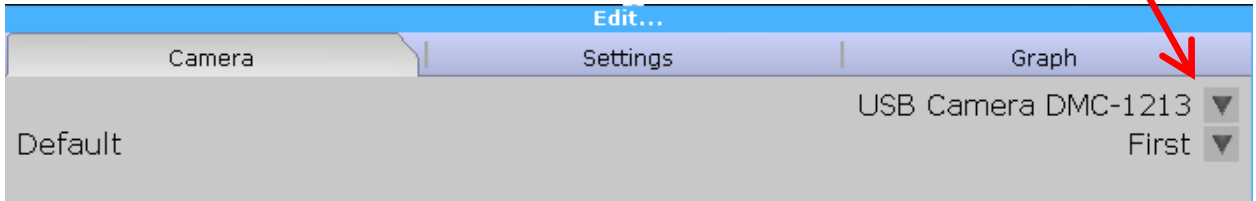

4. Within the second drop-down list, choose **First** to identify this as your primary camera.
5. Click **Apply** and **Close** the edit menu.
6. Now select Camera 2, and click **Edit** to access the edit menu once again.



7. Identify your camera by device name within the first drop-down list and select that camera. If you are only using one camera, select the same camera as before.

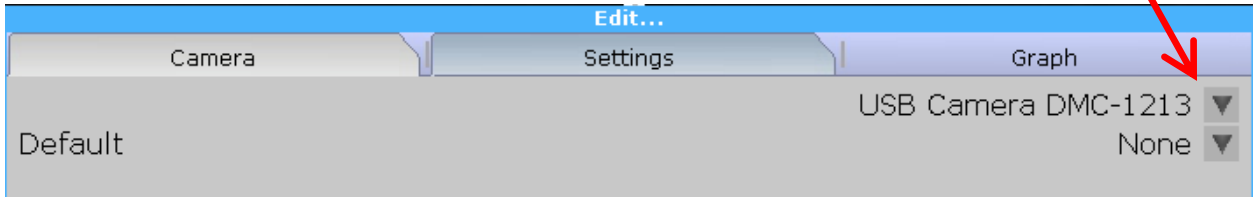

8. Within the second drop-down list, choose **Second** to identify this as your secondary camera. If you are only using one camera, choose **None** instead.
9. Click **Apply** and **Close** the edit menu.
10. Finally, navigate to the **File** Tab and click **Save Workspace** to save these settings.

